

Data Capture revision history

Version 6.0.13 6/17/2024

Minor improvements. Incorrect measurement units for humidity sensors displayed in sensor report.

Version 6.0.12 3/4/2024

Added ability to resend the Fail to Communicate email after 30, 60, 90, 120 minutes if the unit continues to fail to communicate with Data Capture.

Version 6.0.11 8/8/2023

Minor improvement. Data Capture will now send an email every two hours to alert the user if Data Capture fails to communicate with a unit being monitored. Previous versions only sent one email.

Version 6.0.10 4/17/2023

Corrected issue with version .9 when adding the name and comments to the Hi-Low Report.

Version 6.0.9 4/10/2023

Added two fields to allow user to enter their name and comments to the Hi-Low report.

Version 6.0.8 10/27/2022

The sending of an alert email whenever the Data Capture Windows service starts or stops can now be enabled/disabled.

Corrected an issue with the monthly sensor report.

Version 6.0.7 6/20/2022

Corrected error when creating Monthly Sensor report. Error occurred when switching between printers.

Version 6.0.6 6/13/2022

Added comment field to reports

Version 6.0.5

Added option to send status email every 2 hours as well as every 24 hrs.

Increased number of columns to Raw Sensor Data report to save paper.

Version 6.0.4

Updated SMTP security options

Version 6.0.3 10/18/2021

Updated Raw Sensor Data report to include the name of the sensor in the header. Also changed the font color of the data if out of limits; green if below the lower limit, red if above the upper limit.

Version 6.0.2 9/22/2021

Added ability to program eight cell phone numbers for CM705E rev 19 and above.

Version 6.0.1 4/1/2021

Added function to update the time/date chip on all programmed units. This is great for updating for Day Light Savings.

Version 6.0.1 3/10/2021

The big change is writing sensor data to its own individual database ie. tblSensorDataID_XXX.fic. This was done to speed up reports.

Added ability to disable email reminder

Version 5.4.10

Added query return all sensor data for each unit to add these readings to the daily status email.

Data Capture revision history

Added ability to add contact name for each phone number. This is a feature requested by a customer.

Version 5.4.10 1/30/2020


Added support for M408/M409 CO2 monitor.

Version 5.4.10 11/3/2020

Corrected code for downloading log for model VM608E.

Version 5.4.10 10/21/2020

Added code to show IP address in Temperature Guard Units tree.

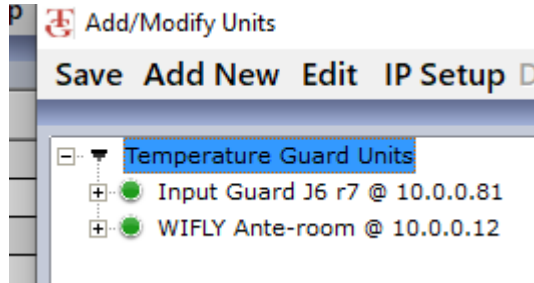

Version 5.4.9 9/22/2020

Corrected the units displayed in the Hi-Lo report for humidity sensor.
Increased font size on all menus.

Version 5.4.8 4/28/20

Added button to check and repair and re-index all database files.

Version 5.4.8 1/2020

Added button to delete unsent alarm emails to the configuration page.
Moved Save, log file buttons to menu bar.

Rev 5.4.7 Compiled and RELEASED 11/8/2019

A local hospital customer requested the ability to edit all unit email addresses at once.
Added menu item under Edit

