

## Data Capture revision history

### **Version 6.0.14 9/27/2024**

Minor improvements. Incorrect measurement units for humidity sensors displayed in sensor report.  
Reduce size of database to increase speed.

### **Version 6.0.13 6/17/2024**

Minor improvements. Incorrect measurement units for humidity sensors displayed in sensor report.

### **Version 6.0.12 3/4/2024**

Added ability to resend the Fail to Communicate email after 30, 60, 90, 120 minutes if the unit continues to fail to communicate with Data Capture.

### **Version 6.0.11 8/8/2023**

Minor improvement. Data Capture will now send an email every two hours to alert the user if Data Capture fails to communicate with a unit being monitored. Previous versions only sent one email.

### **Version 6.0.10 4/17/2023**

Corrected issue with version .9 when adding the name and comments to the Hi-Low Report.

### **Version 6.0.9 4/10/2023**

Added two fields to allow user to enter their name and comments to the Hi-Low report.

### **Version 6.0.8 10/27/2022**

The sending of an alert email whenever the Data Capture Windows service starts or stops can now be enabled/disabled.

Corrected an issue with the monthly sensor report.

### **Version 6.0.7 6/20/2022**

Corrected error when creating Monthly Sensor report. Error occurred when switching between printers.

### **Version 6.0.6 6/13/2022**

Added comment field to reports

### **Version 6.0.5**

Added option to send status email every 2 hours as well as every 24 hrs.  
Increased number of columns to Raw Sensor Data report to save paper.

### **Version 6.0.4**

Updated SMTP security options

### **Version 6.0.3 10/18/2021**

Updated Raw Sensor Data report to include the name of the sensor in the header. Also changed the font color of the data if out of limits; green if below the lower limit, red if above the upper limit.

### **Version 6.0.2 9/22/2021**

Added ability to program eight cell phone numbers for CM705E rev 19 and above.

### **Version 6.0.1 4/1/2021**

Added function to update the time/date chip on all programmed units. This is great for updating for Day Light Savings.

### **Version 6.0.1 3/10/2021**

The big change is writing sensor data to its own individual database ie. tblSensorDataID\_XXX.fic. This was done to speed up reports.

Added ability to disable email reminder

## Data Capture revision history

### Version 5.4.10

Added query return all sensor data for each unit to add these reading to the daily status email.

Added ability to add contact name for each phone number. This is a feature requested by a customer.

### Version 5.4.10 1/30/2020

Added support for M408/M409 CO2 monitor.

### Version 5.4.10 11/3/2020

Corrected code for downloading log for model VM608E.

### Version 5.4.10 10/21/2020

Added code to show IP address in Temperature Guard Units tree.



### Version 5.4.9 9/22/2020

Corrected the units displayed in the Hi-Lo report for humidity sensor.

Increased font size on all menus.

### Version 5.4.8 4/28/20

Added button to check and repair and re-index all database files.

### Version 5.4.8 1/2020

Added button to delete unsend alarm emails to the configuration page.


Moved Save, log file buttons to menu bar.

### Rev 5.4.7 Compiled and RELEASED 11/8/2019


A local hospital customer requested the ability to edit all unit email addresses at once.

Added menu item under Edit

## Data Capture revision history

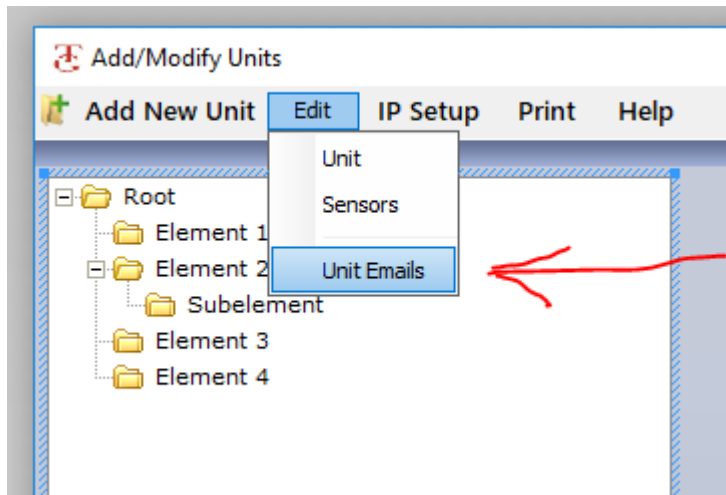

Brings up below screen

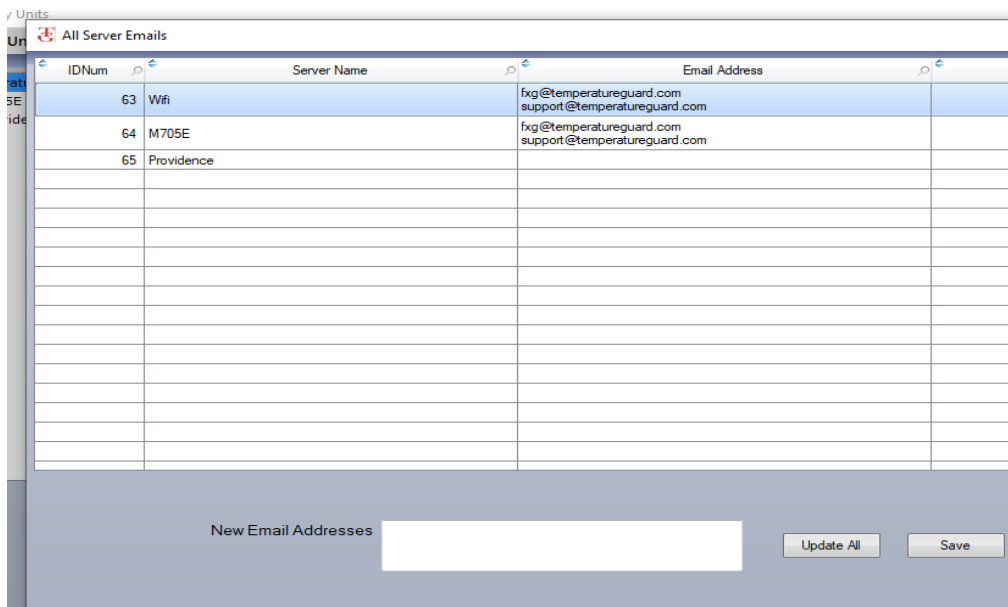

IDNum	Server Name	Email Address
63	Wifi	fxg@temperatureguard.com support@temperatureguard.com
64	M705E	fxg@temperatureguard.com support@temperatureguard.com
65	Providence	

New Email Addresses

Update All Save

### Rev 5.4.6 10/28/2019

Customer requested the Alarm Email Reminder options include 5 minutes.  
The available reminder times are changed depending on the update rate.