

## Data Capture revision history

### Version 6.0.3 10/18/2021

Updated Raw Sensor Data report to include the name of the sensor in the header. Also changed the font color of the data if out of limits; green if below the lower limit, red if above the upper limit.

### Version 6.0.2 9/22/2021

Added ability to program eight cell phone numbers for CM705E rev 19 and above.

### Version 6.0.1 4/1/2021

Added function to update the time/date chip on all programmed units. This is great for updating for Day Light Savings.

### Version 6.0.1 3/10/2021

The big change is writing sensor data to it's own individual database ie. tblSensorDataID\_XXX.fic. This was done to speed up reports.

Added ability to disable email reminder

### Version 5.4.10

Added query return all sensor data for each unit to add these reading to the daily status email.

Added ability to add contact name for each phone number. This is a feature requested by a customer.

### Version 5.4.10 1/30/2020

Added support for M408/M409 CO2 monitor.

### Version 5.4.10 11/3/2020

Corrected code for downloading log for model VM608E.

### Version 5.4.10 10/21/2020

Added code to show IP address in Temperature Guard Units tree.



### Version 5.4.9 9/22/2020

Corrected the units displayed in the Hi-Lo report for humidity sensor.  
Increased font size on all menus.

### Version 5.4.8 4/28/20

Added button to check and repair and re-index all database files.


### Version 5.4.8 1/2020

Added button to delete unsent alarm emails to the configuration page.  
Moved Save, log file buttons to menu bar.

## Data Capture revision history

### Rev 5.4.7 Compiled and RELEASED 11/8/2019

A local hospital customer requested the ability to edit all unit email addresses at once. Added menu item under Edit

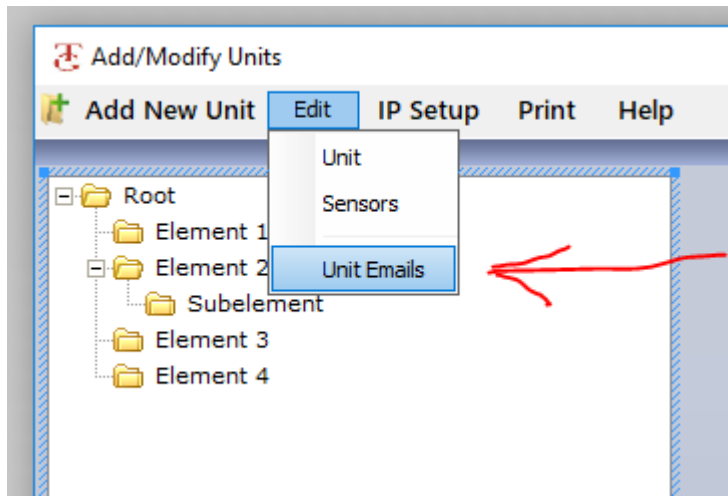

Brings up below screen

The screenshot shows the 'All Server Emails' window with a table of server email addresses. The table has columns for IDNum, Server Name, and Email Address. Below the table is a text input field for 'New Email Addresses' and two buttons: 'Update All' and 'Save'.

IDNum	Server Name	Email Address
63	Wifi	fxg@temperatureguard.com support@temperatureguard.com
64	M705E	fxg@temperatureguard.com support@temperatureguard.com
65	Providence	

### Rev 5.4.6 10/28/2019

Customer requested the Alarm Email Reminder options include 5 minutes. The available reminder times are changed depending on the update rate.

## **Data Capture revision history**